

HOW TO USE LA PIAS EAR TAGS

To prevent infection to the animal's ear decontaminate La Pias Ear Tags and the Ear Tag Applicator using either 70% alcohol or UV irradiation.

1. For the animal's protection it should be anesthetized during this procedure
2. Detach the female portion of the tag with the Applicator (as shown in figure 1).
3. Place this portion into the slot (the red side) of the Applicator (see figure 2 below).



Figure 1

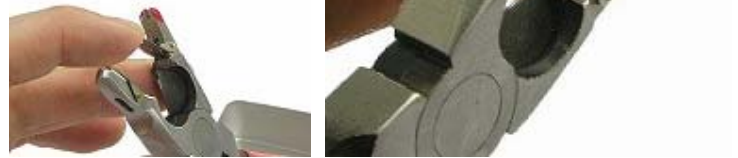


Figure 2

NOTE 1: Make sure that the female portion of the tag is properly placed into the slot before attaching the male portion of the tag on the opposite side of the applicator.

4. Holding the male portion of the tag with forceps place it into the appropriate slot of the Applicator (see figure 3 & 4 below)



Figure 3



Figure 4

Your Distributor is:
Braintree Scientific, Inc.
PO Box 850498, Braintree, MA 02185
781-917-9526
Email: Info@braintreesci.com
Web: www.braintreesci.com

5. Applying force to the handles of the Applicator position the tag in the middle part of the ear as shown in the Figure 5 below. Avoid applying the tag to the ring of cartilage on the ear's edges. Additionally, to avoid bleeding, position the tag away from blood vessels.

NOTE 2: Following application, be sure to check that the tag point has come through the hole properly.



Figure 5



Figure 6 Ear Tag Applicator

Your Distributor is:
Braintree Scientific, Inc.
PO Box 850498, Braintree, MA 02185
781-917-9526
Email: Info@braintreesci.com
Web: www.braintreesci.com