KM CORDLESS
2 Speed Professional Clipper
OPERATION MANUAL

Your Distributor is:
Braintree Scientific, Inc.
PO Box 850498, Braintree, MA 02185
781-917-9526
Email: Info@braintreesci.com
Web: www.braintreesci.com

Can No 94581
©2018 Wahl Clipper Corp.

IMPORTANT SAFETY INSTRUCTIONS

When using an electrical appliance, basic precautions are required to always be followed, including the following: Read all instructions before using this clipper.

DANGER: To reduce the risk of death or injury by electric shock:

1. Do not reach for an appliance that has fallen into water. Unplug immediately from receptacle.
2. Do not use while bathing or in a shower.
3. Do not place or store appliance where it is able to fall or be pulled into a tub or sink. Do not place in or drop into water or other liquid.
4. Always unplug the unit from the electrical outlet immediately after use or after charging the battery.
5. Unplug this appliance before cleaning.

WARNING: To reduce the risk of burns, fire, electric shock, or injury to persons:

1. Do not crush, disassemble, heat above 100°C (212°F), or incinerate the appliance, due to risk of fire, explosion, or burns.
2. Keep this appliance dry.
3. This appliance can be used by children aged from 14 years and above and persons with reduced physical, sensory or mental capabilities or lack of experience and knowledge if they have been given supervision or instruction, by a person responsible for their safety, concerning the use of the appliance in a safe way and understand the hazards involved. Children should be supervised to ensure that they do not play with the appliance. Cleaning and user maintenance shall not be made by children without supervision.
4. The appliance is only to be used with the power supply unit provided with the appliance.
5. Use this appliance only for its intended use as describe in this manual. Do not use attachments not specified by the manufacturer.
6. Never operate this appliance when it has a damaged cord or plug, when it is not working properly, after it has been dropped or damaged, or after it has been dropped into water. Return the appliance to a service center for examination and repair.
7. Keep the cord away from heated surfaces.
8. Never operate this appliance with the air openings blocked or while on a soft surface, such as a bed or couch, where it is possible for the air openings to be blocked. Keep the air openings free of lint, hair, and similar objects.
9. Never drop or insert any object into any opening.
10. Do not use outdoors or operate where aerosol (spray) products are being used or where oxygen is being administered.
11. Do not use this appliance with a blade or attachment that has been damaged or broken, as it has the potential to result in injury.
12. Always attach plug to appliance first, then to outlet. To disconnect, turn all controls to off position (e.g. "0") then remove plug from outlet.
13. During use, do not place or leave the appliance where it is expected to be subject to damage by an animal, or exposed to weather.
14. Cutting blades may become hot after prolonged use.


SAVE THESE INSTRUCTIONS







THIS CLIPPER WAS DESIGNED FOR COMMERCIAL USE ONLY Only use hair clippers to cut human hair. Only use animal clippers to trim animal hair and fur. This appliance is intended for trimming purposes only.

USA and CANADA Lithium Ion Battery Replacement, Repair, and Recycling:

- For battery replacement (US only), send to: Wahl Annex 3001 North Locust Street PO Box 578 Sterling, IL 61081
- Do not attempt to remove the battery.
- Lithium Ion batteries can explode, catch fire, and/or cause burns if disassembled, damaged, or exposed to water or high temperatures.
- If you choose not to send the appliance in for battery replacement, the entire, intact clipper unit must be disposed. Do not attempt to remove the battery for disposal.
- Lithium Ion batteries are subject to local disposal regulations. Many locations prohibit disposal in standard waste receptacles. Check your local laws and regulations.
- For recycling, contact a government recycling agency, waste disposal service, retailer, or visit online recycling sources, such as www.batteryrecycling.com or www.rbc.org.

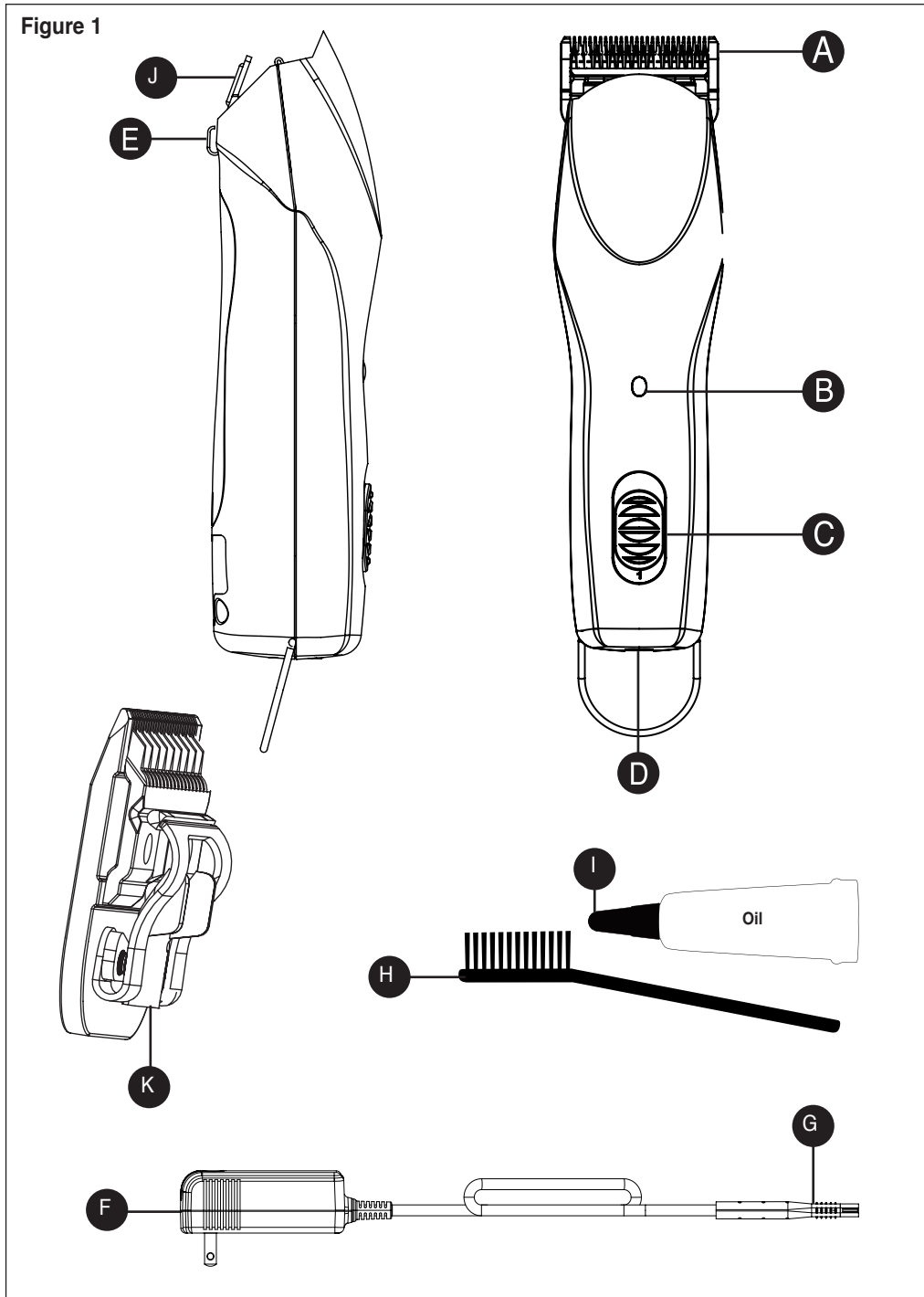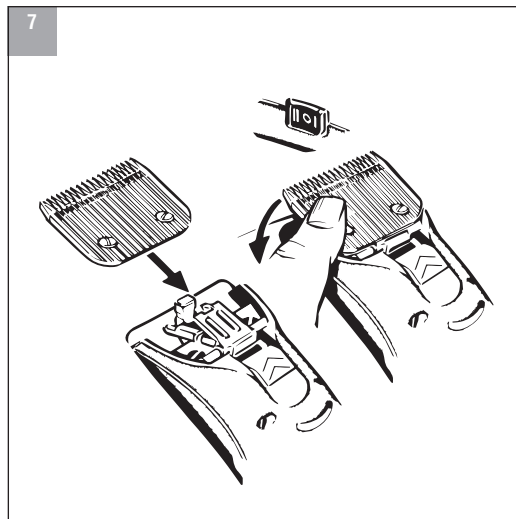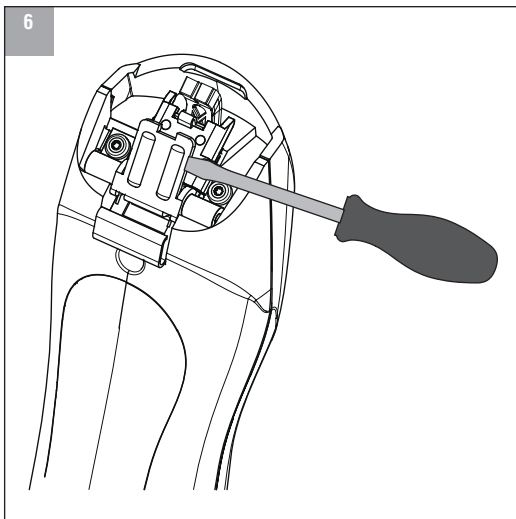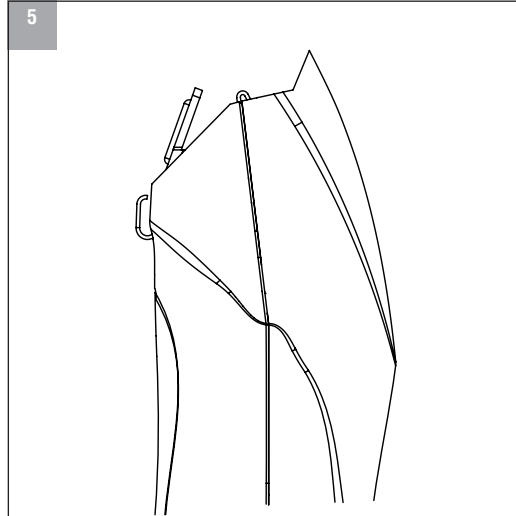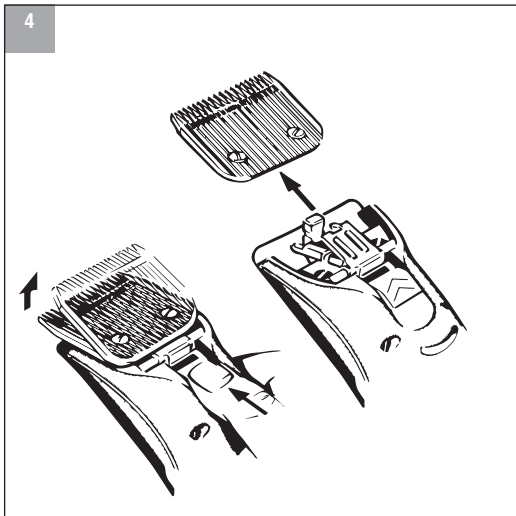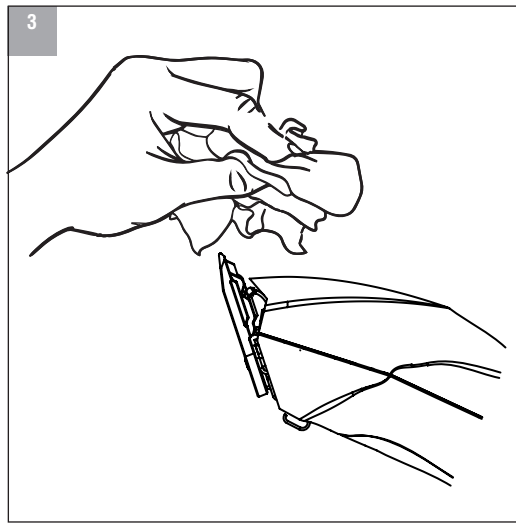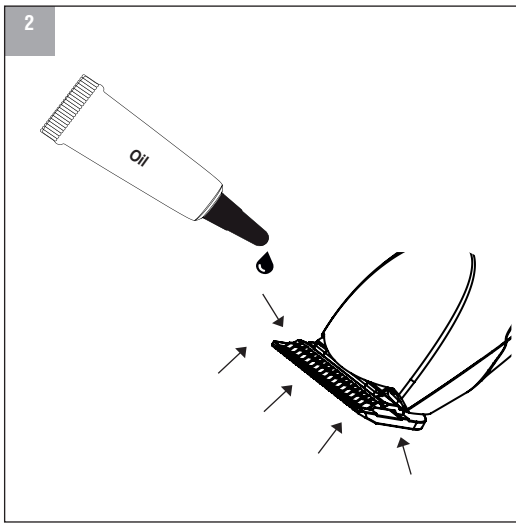Figure 1

DESCRIPTION OF PARTS (FIGURE 1)

- A. Blade set
- B. LED charging indicator
- C. On/off switch
- D. Appliance socket
- E. Blade release button
- F. Appliance plug
- G. Plug-in transformer
- H. Cleaning brush
- I. Blade oil
- J. Blade latch
- K. Blade retainer

DIRECTIONS:

- Charge the clipper battery fully before using for the first time.
- This clipper is equipped with two different operating speeds. Turn the clipper on by sliding the on/off switch upward. Position “1” runs at 3,000 strokes per minute (SPM). Position “2” runs at 3,700 strokes per minute (SPM). After use, switch the clipper off by sliding the on/off switch back down to Position “0”

CHARGING THE BATTERY

- During cordless operation, the LED light will light up solid green. When the green LED light begins to flash, this indicates battery charge is low and the unit will need to be recharged.
- When the battery has fully run down, the clipper will switch itself off automatically.
- To charge, connect the cord to the charging socket on the clipper and then plug the plug-in transformer into an electrical socket.
- When the battery begins to charge, an LED light will blink green to indicate the unit is charging. When the LED light has stopped blinking and remains a constant, unblinking green, the clipper is fully charged and ready for use.
- Full charge time will take approximately 60 minutes. A full charge allows for approximately 120 minutes of clipper run time.
- In the case of low battery, the clipper can be used as a corded unit. However the battery will not begin to charge until the clipper is turned off.
- This clipper includes a “Quick Charge” feature; For example, 10 minutes of charge time will provide 15 minutes of run time.

BATTERY TROUBLESHOOTING

- **Flashing Red LED Light:** The over-current protection circuit has been activated, possibly due to using a dirty or scored blade. The unit will automatically shut down, but can be reset by turning the switch to the off position and then turning it back on. If the problem is not resolved, the unit may stop again right away.
- **Solid Red LED Light:** The battery pack temperature is too high. This could happen if the unit is used under very heavy clipping loads in extreme temperature conditions or while charging the unit in an extreme temperature environment. The unit will be temporarily disabled until the battery temperature returns to a normal level. Do not attempt to recharge the unit until the solid red LED light has turned off. **If a problem persists, please contact Wahl Customer Service at 1-800-PRO-WAHL.**

CLEANING AND CARE

- Do not immerse the clipper in water!
- After each use, use the cleaning brush to remove loose hair from the clipper housing and the blade set.
- The blade set should be oiled frequently (daily) in order to maintain its good cutting performance and thus ensuring a long service life. **Only use Wahl Clipper Blade Oil.** Wahl Clipper blade oil has been specifically formulated for Wahl Clipper blades. Using any other lubricant may cause damage to your blades and clipper and would void your clipper’s warranty.
- Properly oil your blade using the 5 Point Oiling Method (Figure 2). Place 3 drops of oil along the top of the blade teeth at the left corner, right corner, and center of the blade. Turn the clipper to its side and place a drop of oil at the heel of the top blade. Repeat on the other side. Let the clipper run for 30 seconds to allow the oil to work its way into the blade set.
- Wipe away any excess oil. Only wipe the clipper with a soft cloth, which may be dampened slightly. Do not use solvents or abrasives! (Figure 3)
- The blade set may be cleaned with **Wahl Clipper Clini-Clip hygienic spray**. **Do not use any other cleaning agent.** Wahl Clipper Clini-Clip has been specifically formulated for Wahl Clipper blades. Using any other cleaning agent may cause damage to your blades and clipper and would void your clipper’s warranty.
- **Insufficient maintenance of the blade set can reduce the operating time considerably.**
- Care should be used to keep the cord untwisted and undamaged. When storing the clipper, coil the cord. Wrapping the cord around the clipper for storage can cause strain to the cord, which over time, can lead to the loosening of the wires and/or electrical shortages.

CHANGING THE BLADE SET

- Switch off the clipper using the on/off switch.
- Locate the **blade release button** on the backside of the clipper. Press in the blade release button and gently pull the blade away from the clipper. Slide the blade off the **blade latch** (Figure 4).
- To place a new blade on the clipper, ensure that the **blade latch** is in the correct, upright position, not lying flat against the clipper. (Figure 5)
- If the blade latch is not upright, you will need a flathead screwdriver to pry it up. (Figure 6). Use the flathead screwdriver to pry open the blade latch from the side. Do not attempt to pry the blade latch open from the top as there is a chance the drive tip could be damaged.
- Insert the blade onto the blade latch (Figure 7) by sliding the **blade retainer** over the **blade latch**. Before snapping the blade into place, turn the clipper back on. It is important the clipper be running as this ensures the blade will seat properly. Use the palm of your hand to guide the blade to the clipper. There will be an audible “click” when the blade is set in place.

LIMITED 5 (FIVE) YEAR WARRANTY

If your Wahl Product fails to operate satisfactorily within five years from the date of original purchase or receipt as a gift, Wahl will at our option repair or replace it and return it to you FREE OF CHARGE. Our warranty does not cover blades, brushes, or other wear out parts. Please contact Wahl Clipper Corporation's Service and Repair Center located in Sterling, IL at 1-800-PRO-WAHL to receive pre-paid mailing instructions.

Products out of warranty will be repaired at our standard repair charges. This warranty is void and no repair or replacement will be made under this warranty or otherwise if:

1. The product is modified in any manner.
2. The product is repaired by anyone other than Wahl Clipper Corporation or a Wahl Clipper Corporation Authorized Service Center.
3. The product has been subject to unreasonable use.

This warranty gives you specific legal rights and you may also have other rights which may vary from state to state. No other written express warranty is given. Any implied warranty, including any warranty of merchantability, which may arise from purchase or use, is limited to the two-year period in this express warranty. Some states may not allow such limitation, so it may not apply to you. No responsibility is assumed for incidental or consequential damages of any type, or the use of unauthorized attachments. Some states may not allow the exclusion or limitation of incidental or consequential damages, so the limitation or exclusion may not apply to you. Warranty only valid in the United States.

Wahl Clipper Corporation 3001 North Locust Street Sterling, Illinois 61081 • Phone: 1-800-PRO-WAHL Fax: (815) 625-0091

LIMITED 5 (FIVE) YEAR WARRANTY

If your Wahl Product fails to operate satisfactorily within five years from the date of original purchase or receipt as a gift, Wahl will at our option repair or replace it and return it to you FREE OF CHARGE. Our warranty does not cover blades, brushes, or other wear out parts. Please contact Wahl Clipper Corporation's Service and Repair Center located in Sterling, IL at 1-800-PRO-WAHL to receive pre-paid mailing instructions.

Products out of warranty will be repaired at our standard repair charges. This warranty is void and no repair or replacement will be made under this warranty or otherwise if:

1. The product is modified in any manner.
2. The product is repaired by anyone other than Wahl Clipper Corporation or a Wahl Clipper Corporation Authorized Service Center.
3. The product has been subject to unreasonable use.

This warranty gives you specific legal rights and you may also have other rights which may vary from state to state. No other written express warranty is given. Any implied warranty, including any warranty of merchantability, which may arise from purchase or use, is limited to the five-year period in this express warranty. Some states may not allow such limitation, so it may not apply to you. No responsibility is assumed for incidental or consequential damages of any type, or the use of unauthorized attachments. Some states may not allow the exclusion or limitation of incidental or consequential damages, so the limitation or exclusion may not apply to you. Warranty only valid in the United States.

Wahl Clipper Corporation 3001 North Locust Street Sterling, Illinois 61081 • Phone: 1-800-PRO-WAHL Fax: (815) 625-0091

Your Distributor is:
Braintree Scientific, Inc.
PO Box 850498, Braintree, MA 02185
781-917-9526
Email: Info@braintreesci.com
Web: www.braintreesci.com

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