© 2015 Wahl Clipper Corporation
part no. 94677-3402

PROFESSIONAL RECHARGEABLE TRIMMER

IMPORTANT SAFEGUARDS

When using an electrical appliance, basic precautions should always be followed, including the points listed below. Read all instructions and safeguards before using.

DANGER

To reduce the risk of injury or death by electric shock:

1. Do not reach for an appliance that has fallen into water. Unplug immediately from the outlet.
2. Do not use while bathing or in a shower.
3. Do not place or store appliance where it can fall or be pulled into a tub or sink. Do not place in or drop into water or other liquid.
4. Except when charging, always unplug this appliance from the electrical outlet immediately after using.
5. Always unplug this appliance before cleaning.

WARNING

To reduce the risk of burns, fire, electric shock or injury to persons:

- This appliance can be used by children aged from 14 years and above and persons with reduced physical, sensory or mental capabilities or lack of experience and knowledge if they have been given supervision or instruction concerning use of the appliance in a safe way and understand the hazards involved. Children shall not play with the appliance. Cleaning and user maintenance shall not be made by children without supervision.
- Warning: Keep the Product Dry. It is essential to PREVENT electrical equipment from coming into contact with water or other liquids.
- The appliance is only to be used with the power supply unit provided with the appliance.
- Only use hair clippers to cut human hair.
- Only use animal clippers to cut animal hair and animal fur.

Your Distributor is:

Braintree Scientific, Inc.

PO Box 850498, Braintree, MA 02185

781-917-9526

Email: Info@braintreesci.com

Web: www.braintreesci.com

- The appliance is intended for trimming purposes only.
 - **WARNING:** Cutting blades may become hot after prolonged use.
 - Use this appliance only for its intended use as described in this manual.
- Do not use attachments not recommended by the manufacturer.
- The appliance is intended for trimming purposes only.
 - Never operate this appliance if it has a damaged cord or plug, if it is not working properly, if it has been dropped or damaged, or if it has been dropped into water. Return the appliance to Wahl Clipper Corporation for examination and repair.
 - Keep the cord away from heated surfaces.
 - Insert only recharging transformer plug into charger receptacle.
 - Do not use outdoors or operate where aerosol (spray) products are being used or where oxygen is being administered.
 - Do not use this appliance with a damaged or broken comb nor with teeth missing from the blades, as injury may occur.
 - During use, do not place or leave appliance where it may be damaged by an animal or exposed to weather.
 - Always attach plug to appliance first, then to outlet. To disconnect, turn unit "OFF" then remove plug from outlet.
 - Do not crush, disassemble, heat above 100°C (212°F), or incinerate the appliance, due to risk of fire, explosion, or burns.

SAVE THESE INSTRUCTIONS FOR COMMERCIAL USE ONLY

Description of parts

- A Blade set
- B On/off switch
- C Charging indicator
- D Charging stand
- E Blade guard
- F Cleaning brush
- G Oil for blade set
- H 97581 series plug-in transformer (shape may differ, depending on country of origin)

Charging the battery

Before the appliance is used for the first time, it should be charged for 2 hours. It then takes 1 hour to fully recharge the battery again.

Connect the plug-in transformer to the appliance socket, and then to the mains socket.

The appliance can also be charged via the charging stand. Connect the plug-in transformer to the charging stand's socket. Place the appliance in the charging stand so that charging contact is made.

During charging, the charging indicator is illuminated. The charging indicator will begin to blink when almost fully charged and may shut off when fully charged.

Note: If the appliance is fully charged and is placed in the stand to charge, the charging indicator may not turn on. This does not indicate the appliance can no longer charge, it means the appliance does not need to draw electricity because it is already fully charged.

Mains operation

Connect the plug-in transformer to the appliance socket (Fig. 2). Connect the plug-in transformer to the mains socket.

Switch on the appliance using the on/off switch (Fig. 1a) and, after use, switch it off again (Fig. 1b).

Cleaning and care

Oil the blade set (Fig. 3).

Do not immerse the appliance in water!

After each use, use the cleaning brush to remove cut hairs from the housing opening and from the blade set (Fig. 5). Remove the blade set to do this (Fig. 4a).

The charging contacts on the appliance and on the charging stand should always be kept clean. Only wipe the appliance with a soft cloth, which may be dampened slightly. Never use solvents or scouring agents!

The blade set can be cleaned with hygienic spray order no. 4005-7051.

For a long-lasting, good cutting performance, it is important to oil the blade set frequently (Fig. 3). Only use blade set oil order no. 0230-1070 (118 ml) or no. 3310-230 (4oz.). Blade set oil and hygienic spray is available from your dealer or our Service Centre.

If the cutting performance deteriorates after long use despite regular cleaning and oiling, the blade set should be replaced.

Use the blade guard when the appliance is not in use.

Always store the appliance with the blade guard in place.

Changing the blade set

Switch off the appliance using the on/off switch (Fig. 1b).

The blade set can be removed by sliding it away from the housing in the direction of the arrow (Fig. 4a).

To replace the blade set, position it with the hook in the opening of the housing and press it onto the housing until it clicks into position (Fig. 4b).

This marking indicates that this product should not be disposed with other household wastes throughout the EU. To prevent possible harm to the environment or human health from uncontrolled waste disposal, recycle it responsibly to promote the sustainable reuse of material resources. To return your used device, please use the return and collection systems or contact the retailer where the product was purchased. They can take this product

for environmental safe recycling.

Lithium Battery Replacement, Removal and Disposal

Lithium-Ion Battery Replacement

For battery replacement the entire, intact trimmer should be sent to: Wahl Customer Services (the address can be found on the top of this booklet) or taken to a suitably qualified professional. The battery should only be replaced by the original manufacturers spare, for safety reasons a substitute should not be used.

Do not attempt to remove the battery. Lithium-Ion batteries can explode, catch fire, and/or cause burns if disassembled, damaged, or exposed to water or high temperatures.

Lithium Battery Disposal

Do not dispose of this product or the Lithium battery in the normal household waste. The product should be returned to Wahl Customer Services where we offer a take back service or taken to a local collection facility where they will be dealt with safely and sent for the appropriate recycling.

SERVICE AND REPAIR

-This appliance contains no user-serviceable parts.

-Any servicing should be done by Wahl Service and Repair or by an Authorized Service Center.

-Do not return this product to the store where purchased. If you have problems with your unit, send it prepaid with a note explaining the problem to:

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