

Super Pocket Pro OPERATING INSTRUCTIONS

BATTERY INSTALLATION/REPLACEMENT

1. Remove battery compartment cover from bottom of trimmer by sliding it off.
2. Insert one AA alkaline battery into trimmer as shown.
3. Replace battery compartment cover to its original position.

NOTE: To maximize battery life, trimmer should be shut "OFF" immediately after each use.

HOW TO USE THE TRIMMER ATTACHMENTS

Your WAHL Pocket Pro trimmer includes two trimming attachments - medium and long cut. To use either one of these attachments, simply insert one side of attachment over the front of the blades and then push the other side back until both sides are locked into position. To remove attachment, pull tab away from one side to release attachment from trimmer. Use somewhat of a rolling pressure action, with your thumb on top and index finger underneath.

CAUTION: Make sure that the guide is securely in place on blades before operating unit. The two small plastic tabs inside the guide should be OVER the blade and locked in the blade groove.

This trimmer offers an easy way to trim around eyes, ears and legs without shaving hair completely off. To trim the length you want, just choose the attachment and cutting directions indicated in the chart below. After selecting the length and trimmer guide desired, begin trimming your animal.

Desired Length	Attachment	Direction of Cut
.8mm (1/32")	None	With lay of hair
1.6mm (1/16")	Short	Against lay of hair
3mm (1/8")	Short	With lay of hair
5mm (3/16")	Long	Against lay of hair
6mm (1/4")	Long	Against lay of hair

WHERE TO BEGIN:

Begin by bathing your animal. Dirt or sand particles in an animal's coat will quickly dull trimmer blades. Rinse well and allow the coat to dry completely. Comb through the coat to remove any tangles and make trimming as easy as possible. This unit is intended for trimming and touch-ups. For complete clipping of body hair, you will need a more powerful clipper.

SMALL ANIMALS:

A little trimming will really enhance a dog's appearance. Carefully begin trimming the legs and feet according to the style desired. The legs may be trimmed while your animal is sitting on a table, with the leg in your hand or while the animal is standing. Trimming with the grain or lay of hair will give a slightly longer cut. Be careful when trimming around joints and projections. Continue by trimming around the muzzle, chin and ears. Hold your dog's muzzle with your free hand and trim around the face and chin. To trim the ears, lay the ears in the palm of your hand. Trim from the top of the ears downward, clipping the outside of the ear first and then the inside. Hold the ear with your free hand and trim the sides.

OILING BLADES:


For best results, put a drop of Wahl Clipper Oil on blades periodically, or approximately once a month. Motor bearings are permanently lubricated. Do not attempt to oil.

NOTICE-DO NOT OVER OIL

This trimmer requires very little oiling. When necessary to oil, turn trimmer "ON" hold the unit with blades facing DOWN, and squeeze one drop of the oil onto blades. Wipe off excess oil with soft cloth.

T-Blade Installation Instructions

1. Remove the old blade set by removing the two screws.(1)
2. Clean excess hair from open end of trimmer with a small brush.
3. Position cam follower (2) with flat side up into top blade(3) holes.
4. Align each spring hook(4) with the slots in the front on each side of the cam follower(2). Then push the top blade(3) and cam follower(2) assembly back from the heel area of the top blade(3) and hold it back. Make sure that the spring hooks (4) stay inside the shallow slots of the cam follower(2) in the front.
5. While holding the top blade(3) down, position bottom blade(5) over top blade(3) aligning screw holes.
6. Holding the blade set in place, replace the screws and tighten.
7. Turn the trimmer "ON" to check for proper back and forth blade motion.
8. If the blades do not operate properly, loosen screw and slightly lift blades set to check alignment, as discussed in Step 4. Tighten screws and try again.

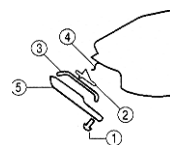

DO NOT ATTEMPT ANY ADJUSTMENTS WHILE TRIMMER IS IN OPERATION. ALWAYS SHUT MOTOR "OFF" FIRST.

ONE-YEAR LIMITED WARRANTY

If your Wahl product fails to operate satisfactorily within one (1) year from the date of original purchase or receipt as gift, Wahl will, at our option, repair or replace it and return it to you FREE OF CHARGE. Our warranty does not cover blades, which are "wear out" parts. Please mail your product prepaid to Wahl Clipper Corporation, Wahl Annex, 3001 N. Locust Street, Sterling Illinois, 61081, Attention: Repair Department, with a note describing the defect and the date of original purchase or receipt as gift (if in warranty). Products out of warranty will be repaired at our standard repair charges. This warranty is void and no repair or replacement will be made under this warranty or otherwise if : 1) The product is modified in any manner. 2) The product is repaired by anyone other than Wahl Clipper Corporation. 3) The product has been subject to unreasonable use.

This warranty gives you specific legal rights and you may also have other rights which may vary from state to state. No other written express warranty is given. Any implied warranty, including any warranty of merchantability which may arise from purchase or use, is limited to the one year period provided in this express warranty. Some states may not allow such limitations, so it may not apply to you. No responsibility is assumed for incidental or consequential damages of any type, or for the use of unauthorized attachments. Some states may not allow the exclusion or limitation of incidental or consequential damage, so the limitation or exclusion may not apply to you. Warranty valid in U.S.A. only. Outside U.S.A., check with store for Wahl authorized distributors' warranty.



Your Distributor is:

Braintree Scientific, Inc.

PO Box 850498, Braintree, MA 02185

781-917-9526

Email: Info@braintreesci.com

Web: www.braintreesci.com