

Dyets Powder Feeder Jar Setup Instructions



- Prevents tipping or moving of Dyets Feeding Jar
- 100% Stainless Steel for Easy Cleaning and Durability
- Versatile Design Allows for Placement in Any Cage Corner
- Set contains jar, lid, follower and holder

Place the holder (base) in the cage under the bedding (no bedding inside the ring)

Fill the jar with powdered food and lay the follower on top of the food.

Slide the lid onto the jar with the bent tabs down and twist to tighten.

Place filled jar in holder (base)

Your Distributor is:
Braintree Scientific, Inc.
PO Box 850498, Braintree, MA 02185
781-917-9526
Email: Info@braintreesci.com
Web: www.braintreesci.com