

Hydrophobic sand that keeps the animal's urine on top, making sample collection easy.

Suitable for most urinalysis tests as a metabolic cage replacement!

LabSand is the name of our patented hydrophobic sand that keeps animal urine on top, making sample collection stress-free. The sand has a natural non-toxic repellent coating. It allows veterinarians, technicians and scientists to passively collect a urine sample without resorting to collection methods like metabolic caging. The kit comes with everything you need to collect a urine sample.

Directions for Use:

Simply spread the contents of a LabSand bag onto a clean surface, and let the animal freely urinate on the sand. The urine will stay beaded on top, allowing you to collect the sample with the provided pipette and vial. Conversely you can also dip a testing strip directly into the droplet while it is on the sand.

LabSand has been validated by several pharmaceutical companies. To view our validation studies please visit www.labsand.com. LabSand is based on the proprietary technology used in Kit4Cat, a veterinary product widely used for feline urine samples in 44 countries worldwide.



Noninvasive



100% Natural



Easy to Use




Suitable for testing all common parameters





Benefits for Scientists, Researchers, & Technicians:

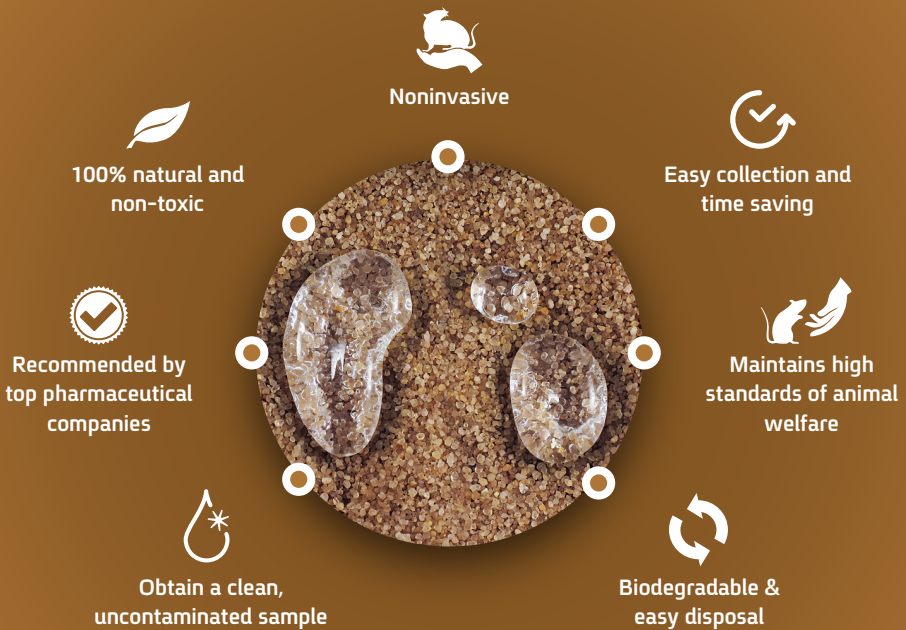

Your Distributor is:
Braintree Scientific, Inc.
PO Box 850498, Braintree, MA 02185
781-917-9526
Email: Info@braintreesci.com
Web: www.braintreesci.com