

Inserting and Changing Burrs Addendum



Your New Micro Drill comes with a plastic insertion/removal burrs tool to help you insert and remove Burrs from the drill's metal chuck



To insert burrs, always place your index finger firmly on the chuck's dome, next insert the burr to the opening and press it in. The tip of the burr should stick out about 20 mm



To remove the burrs, place your index finger firmly on the chuck's dome, insert the removal tool as shown, then pull the burr out. If the metal chuck opening becomes loose-fitting, use a small pair of fine pliers, to squeeze tight the opening. If you over pressed the opening of the metal chuck, then use the other end of the burr to insert and re-open it to the burr's diameter

Your Distributor is:
Braintree Scientific, Inc.
PO Box 850498, Braintree, MA 02185
781-917-9526
Email: Info@braintreesci.com
Web: www.braintreesci.com