

IMPORTANT!!!! PLEASE READ BEFORE USING:

CARE AND USE NOTES

RIS (Rodent Intubation Stand)

Constructed from durable polycarbonate, this platform provides the perfect stage for intubating mice and rats of all sizes. The stand is constructed from a single piece of polycarbonate plastic with a seamless design. The top of the stand provides a hand rest for stabilization during intubation that is designed to support significant weight.

OPERATION

The animal is suspended from the incisor wire on the angled side of the stand. Three mounting positions are available to customize the placement of the animal on the stand in order to meet the user's needs. Two sets of posts are provided with each stand for installation of the incisor wire and the optional head positioning wire. Each post has three positions for the height adjustment to optimally accommodate the user and the animal. Each post is also fitted with a slit for securing the end of the wire making the wire easy to install and easier to remove.

CLEANING

The compact seamless design and its construction from durable polycarbonate make this stand easy to clean and will not crack with the use of harsh cleaning agents as is common with similar stands constructed from acrylic.

STORAGE

The stand is lightweight and portable so it can be stored in the most convenient location possible. The posts are removable and should be stored with the stand.



Your Distributor is:
Braintree Scientific, Inc.
PO Box 850498, Braintree, MA 02185
781-917-9526
Email: Info@braintreesci.com
Web: www.braintreesci.com